

Buhl VP20ADPT Microscope Adapter

Installation Instructions

Thank you for purchasing the VP20ADPT Microscope Adapter for the Buhl VP20M visual presenter. This adapter installs in seconds to your VP20M.



1. Slide the end of the VP20ADPT on to the end of your VP20M's lens as shown:
2. Install the 3 included thumbscrews to secure the VP20ADPT to the VP20M lens
3. Aim VP20M with Microscope Adapter down the eyepiece of any microscope and adjust the focal length on both the microscope and VP20M lens until picture becomes clear.

The VP20M's focus is adjusted by rotating the end of the lens.

For Questions and Support please contact us at support@buhl-ind.com or visit our website at www.buhl-ind.com.